


搅拌液体

连接电源前，确保当地电源电压匹配搅拌器要求。

1. 将搅拌棒放入容器并填充需要搅拌的液体。
2. 将容器放置在工作盘中心位置。
3. 顺时针旋转旋钮打开搅拌器并调节搅拌速度。
4. 如果搅拌棒运行不畅，略微移动容器，确保其始终位于容器与工作盘的中心位置。

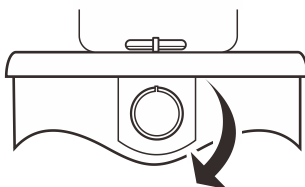

技术参数

最大搅拌容量	2000 mL (H ₂ O)
搅拌速度	0 至 1250 转/分钟
工作面板直径	145 mm (5.7 in.)
工作面板材质	PC
电源类型	AC 220V/50Hz
外形尺寸	185 (Ø) × 75 mm (7.28 × 2.95 in.)
重量	0.6 kg (1.3 lb)

Disposal

This product is required to comply with the European Union's Waste Electrical and Electronic Equipment (WEEE) Directive 2002/96/EC and may not be disposed of in domestic waste. Please dispose of product in accordance with local regulations at the collecting point specified for electrical and electronic equipment.



Warranty

The warranty period for stirrer is one year from the date of shipment. Above warranty does not cover the stir bar. Out of warranty products will be repaired on a charged basis. The warranty on your stirrer shall not apply to defects resulting from:

- Improper or inadequate maintenance by customer
- Unauthorized modification or misuse
- Operation outside of the environment specifications of the products

For more information, please contact the supplier.